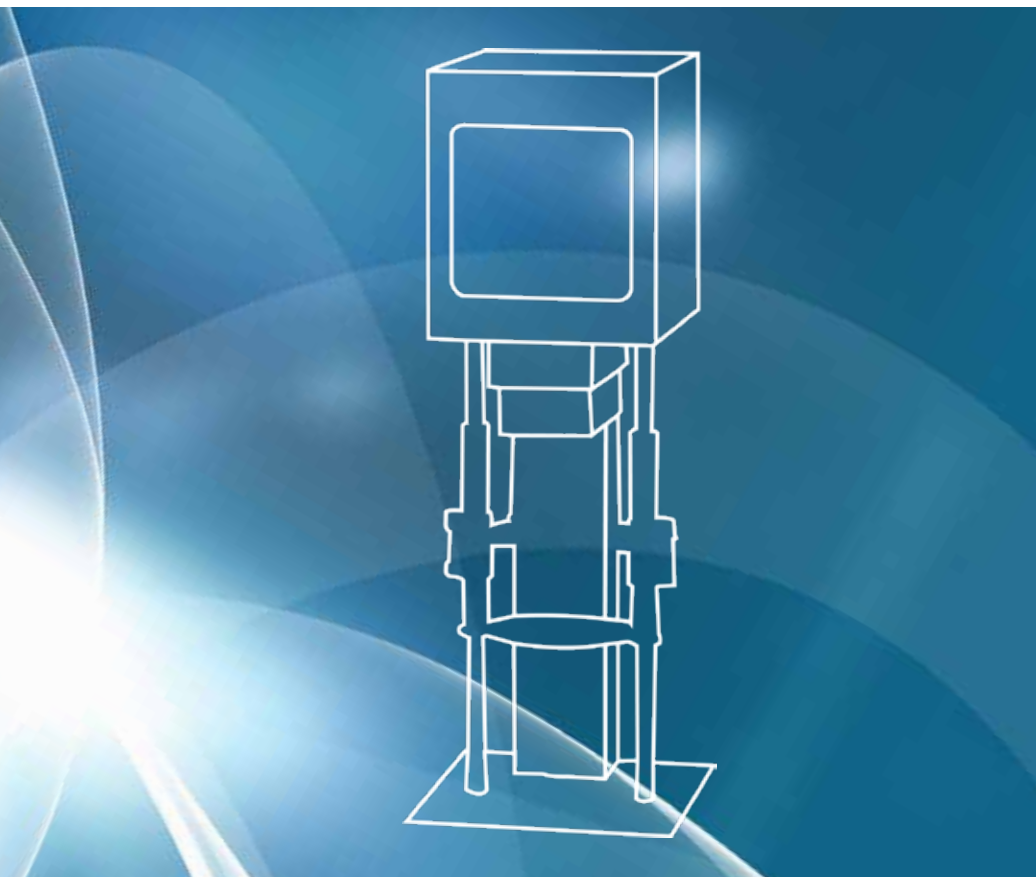
TOPITEC® AUTOMATIC

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im Sinne der

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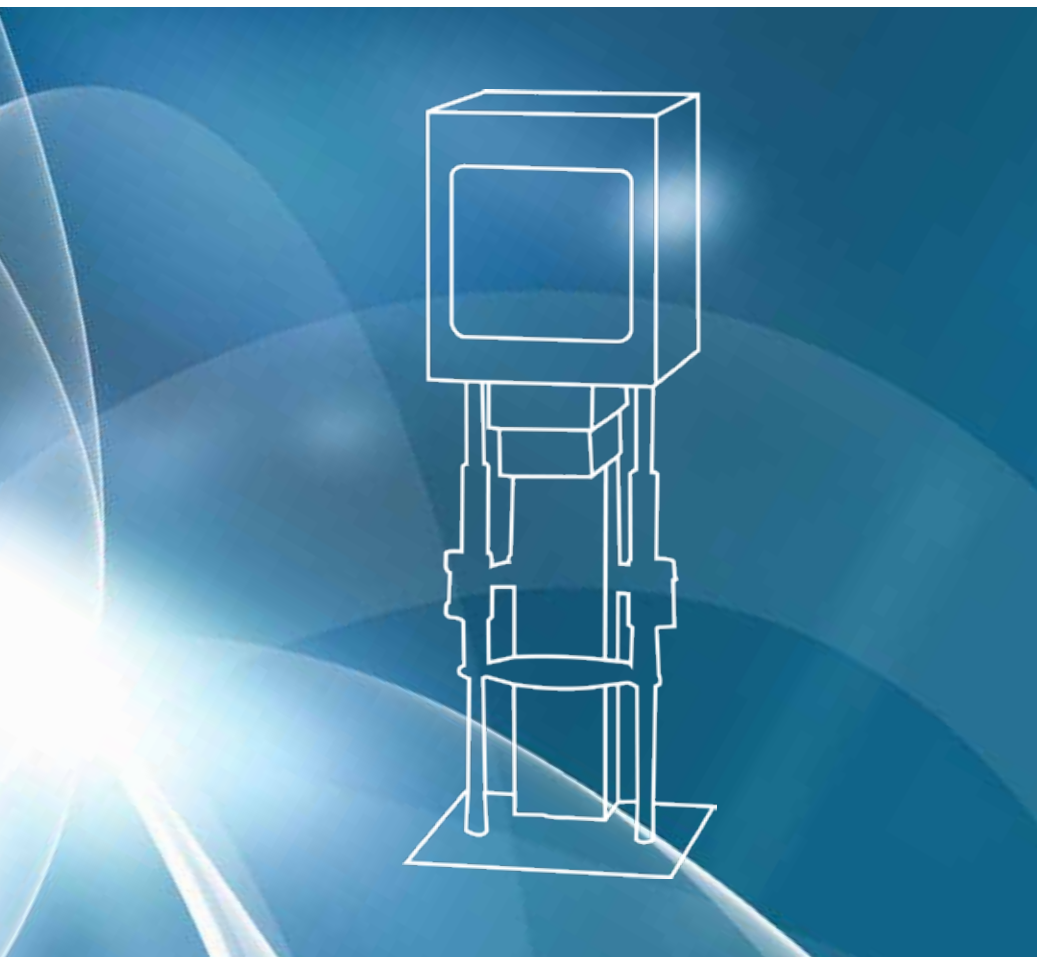
DECLARATION OF CONFORMITY

In accordance with

- **EC Machinery Directive 2006/42/EC**
- **EU Low Voltage Directive 2014/35/EU**
- **EU EMC Directive 2014/30/EU**
- **EU RoHS Directive 2011/65/EU**

We herewith declare that the design of

product: TOPITEC® AUTOMATIC
S/N; machine no. 27XXX-01

as brought onto the market by us conforms to the
DIN EN standards listed below. Harmonised standards:

Directive / standard	Title
2014/30/EU	EC Directive: EMC Directive valid from 01/04/2016
DIN EN 61000-4-2	Testing of immunity to interference from the discharge of static electricity
DIN EN 61000-4-4	Testing of immunity to interference from quick transient electrical disturbance variables/burst
DIN EN 61000-4-5	Testing of immunity to surge voltages
DIN EN 61000-4-6	Immunity to conducted disturbances induced by high-frequency fields
DIN EN 61000-6-3	Emitted interference for residential, commercial and light-industrial environments
2014/35/EU	EC Directive: Low Voltage Directive valid from 01/04/2016
DIN EN 61010-1:2011	Safety requirements for electrical equipment for measurement, control, and laboratory use – Part 1: General requirements
2006/42/EC	EC Directive: Machinery Directive valid from 29/12/2009
DIN EN ISO 12100:2011	Safety of machinery – General principles for design – Risk assessment and risk reduction
DIN EN 349 :2008	Safety of machinery – Minimum gaps to avoid crushing of parts of the human body
2011/65/EU	EC Directive: RoHS valid from 03/01/2013

- This Declaration applies exclusively to the product as delivered.
- The basic requirements as regards health and safety as laid down in Annex I of Machinery Directive 2006/42/EC are fully complied with.
- The special technical documentation according to Annex VII A has been compiled and shall be made available in electronic format to the relevant state authorities on request.

Note

This operating manual forms an integral part of the scope of delivery of the mixing system. A copy of this manual must be accessible at all times to operating and maintenance personnel. All safety instructions in this document must be strictly adhered to. When selling the mixing system, the operating manual must be handed over to the new owner.

You can find lots of useful tips about TOPITEC® at www.topitec.de. The appropriate formulation manual is also available for you to download.

TRANSLATION

Machines shipped to EFTA countries are accompanied by a translation of the operating instructions into the respective language. In the event of discrepancies between the translation and the original (German) version of the manual, the German version is deemed valid. If you have any queries, contact the manufacturer.

COPYRIGHT

The copyright in this document is owned by WEPA Apothekenbedarf GmbH & Co KG. The reproduction or disclosure of the content of this user manual, in any format whatsoever, is only permitted with the prior written consent of WEPA Apothekenbedarf GmbH & Co KG. WEPA Apothekenbedarf GmbH & Co KG shall not be liable for damage caused in connection with the use of this manual.

Reproduction of this document, its use and the communication of its content to a third party are only permitted with the express consent of WEPA Apothekenbedarf GmbH & Co KG. Any breach will result in legal action for damages. All rights reserved.

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TABLE OF CONTENTS

1 Content

1.1 Table of contents	4-6
1.2 Introduction	7
1.3 Service	8

2 Device diagram and proper use 9

2.1 Diagram of TOPITEC® AUTOMATIC	9
2.2. Proper use	10-11
2.3 Technical data	12-13

3 Safety 14

3.1 Safety instructions and notes	14-16
3.2 Built-in safety systems	17-18
3.3 Duties of device owner	19-20
3.4 Safety tests and safety checks	21

4. General safety instructions 22

4.1 Dangers	22
4.2 Danger areas of TOPITEC® AUTOMATIC	22
4.3 Operating and maintenance personnel	23
4.4 Replacement of parts	24
4.5 Shut-down procedure	25

5 Standard equipment 26

5.1 Scope of delivery	26
5.2 Standard equipment	26
5.3 Installation	27
5.4 Transport	28
5.5 Installation and assembly	29
5.6 Commissioning	29

6 Operation	30
6.1 Connection to power supply, interfaces	30-31
6.2 Safety bracket	32
6.3 Control elements of TOPITEC® AUTOMATIC	33
6.4 Operating modes and control	34
7 Pricing	48-49
8 Accessories of TOPITEC® AUTOMATIC	50-53
9 Cleaning / maintenance / inspection / function testing	54
9.1 Maintenance	57
9.2 Cleaning	56-57
9.3 Inspection intervals – function testing	58
9.4 Warning signs attached to TOPITEC® AUTOMATIC	58
9.5 Warranty and liability	59
10 Troubleshooting	60
10.1 Practical tips – instant remedy	60-61
11 Emergency-stop	62
12 Disposal	62
13 Notes	63
14 Index	64-65

TABLE OF CONTENTS

1.2 List of figures

Fig. 1-1 Control elements of TOPITEC® AUTOMATIC	9
Fig. 1-2 Scope of delivery and colours of TOPITEC® AUTOMATIC	26
Fig. 1-3 Interfaces, connection to power supply, fuse holders TOPITEC® AUTOMATIC	30
Fig. 1-4 Safety bracket of TOPITEC® AUTOMATIC	32
Fig. 1-5 aponorm® packaging units for TOPITEC® AUTOMATIC	48-49
Fig. 1-6 Accessories of TOPITEC® AUTOMATIC	50-53

1.3 List of tables

Table 1 Test intervals for safety systems	17
Table 2 Weights of transport unit	28
Table 3 Warning signs attached to TOPITEC® AUTOMATIC	58

INTRODUCTION

The production of preparations by pharmacists for the treatment of dermatological disorders is bound to remain an important element of drug treatment. In this context, the preparation of medicine for external application plays an important role.

The move from manual grinding with mortar and pestle to semi-automated and automated production methods offers many advantages as regards the quality and the safety of pharmaceutical products.

With the innovative TOPITEC® AUTOMATIC, you now have the option to perform many different mixing processes and to store relevant parameters so that they can be called up at any time for other batches. Whether you use the TOPITEC® jar to mix ingredients for a single batch or work with the larger stock preparation container available as an optional accessory: The TOPITEC® AUTOMATIC can save your settings and formulations, if that is useful to you.

To talk to our customer support specialists, call:

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CUSTOMER SERVICE

You can find the quick help you need here.

News and tips

The latest information and many useful tips in connection with the production of medicines with the TOPITEC® mixing system are published on our website at www.topitec.de. Please visit our website and share your experiences with the device with other users!

Galenic queries concerning TOPITEC® equipment

Our qualified pharmaceutical specialists assist you in all matters regarding formulations and applications that might arise in the use of your TOPITEC® mixing system.

Telephone +49(0) 2624 107-145

info@topitec.de

www.topitec.de

Technical support for TOPITEC® equipment

Our technical support staff will be able to provide you with immediate telephone assistance in the event of technical malfunctions on the TOPITEC® mixing systems. In Germany, we operate a technical call-out service.

Please call us on the following number:

Telephone +49(0) 2624 107-361

info@topitec.de

www.topitec.de

CONTROL ELEMENTS

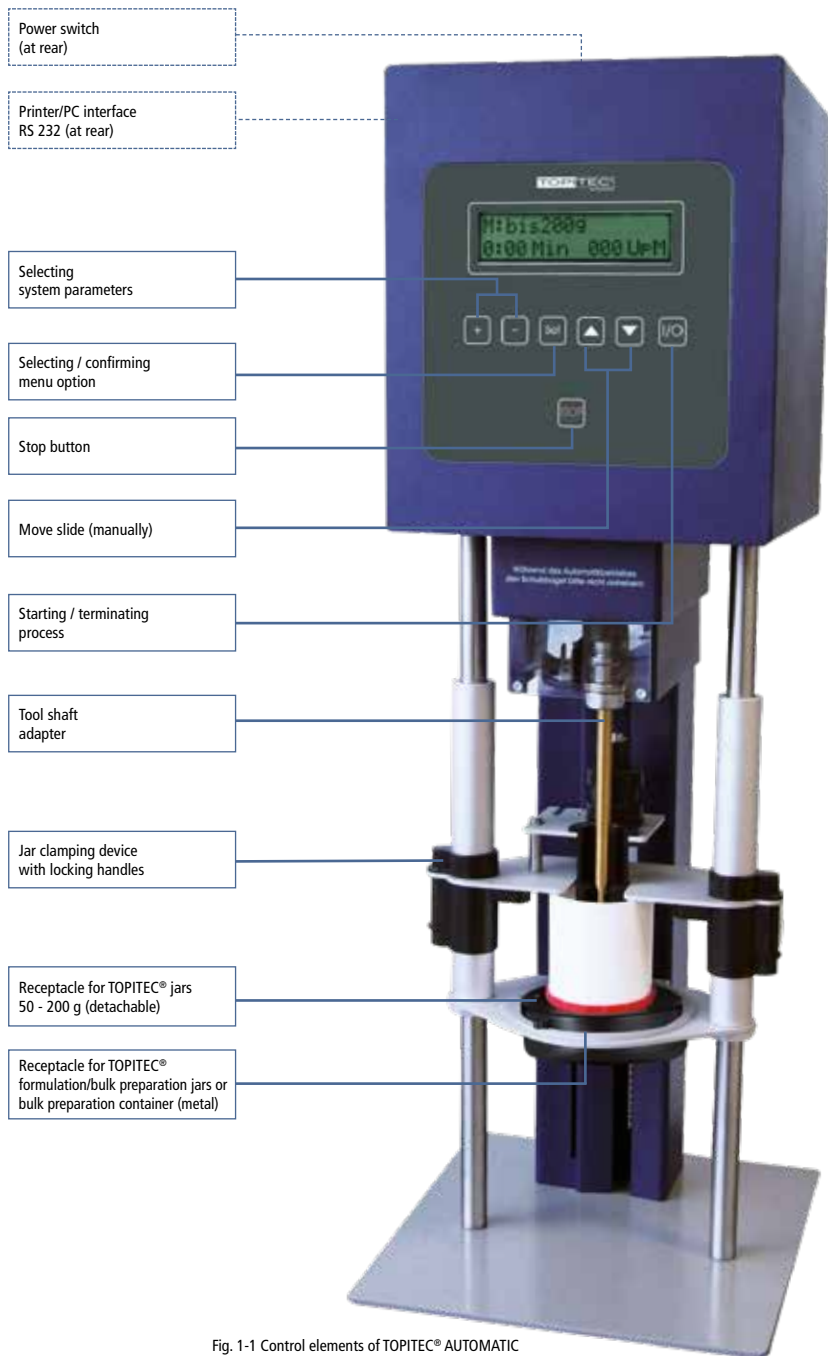
Fig. 1-1 Control elements of TOPITEC® AUTOMATIC

2.1. Proper use / product description

- The TOPITEC® AUTOMATIC is designed for the production of semi-solid formulations (medicines, ointments, creams, cosmetic products) in pharmacies.
During production, the jar containing the ingredients is moved up and down while the product inside the jar (TOPITEC® jar) is stirred by means of the mixer (tool shaft). The preparation of the product (adding ingredients in the correct sequence to the jar) takes place outside the device.
Instruction by WEPA or WEPA sales partner.
- The TOPITEC® AUTOMATIC has been developed and designed exclusively for commercial applications. Any private use of the TOPITEC® AUTOMATIC is forbidden.
- The TOPITEC® AUTOMATIC is not designed for operation in potentially explosive atmospheres (vapour or dust; ATEX; observe information in safety data sheets) or within a designated ATEX zone as defined by the device owner.

2.2 Proper use / application

Mixing of prepared and sealed jars containing pH neutral pastes, ointments, creams or powders that are not radioactive, explosive or flammable. The device is designed for use in pharmacies and operation by suitably instructed personnel. The mixing of ingredients must not result in explosive substances.

An instructed person (operator) is a person with a pharmaceutical qualification who is older than 18 years of age and has been instructed in the use of the device.

The substances, materials and media to be used for processing in the mixing system are to be sourced by the device owner, who is responsible for the proper use of these substances. The device owner is also responsible for the correct handling of these substances, materials and media, with due regard to any associated risks. The device owner must provide safety and disposal instructions for staff. Strictly observe the safety data sheets provided by the manufacturers of the materials / media.

For safety and quality reasons, the device must only be used in conjunction with original TOPITEC®/aponorm® containers. These containers are not gas-tight, so substances might leak out. The device owner must ensure that all personnel wear the required personal protective equipment. Compliance must be monitored.

Operating personnel must comply with all statutory regulations and observe the specific requirements for the processed substances.

Information regarding proper hygiene procedures can be found in the ApBetrO, the hygiene guidelines published by Gesellschaft für Dermopharmazie GD for the preparation of sterile pharmaceutical products, the BAK Guidelines, statutory regulations, etc.

Warning

The mixing system must only be used for the purposes listed above. Any other use is deemed improper and is thus prohibited. The mixing system must not be modified in any way, unless written consent has been obtained from the manufacturer.

The manufacturer shall not be liable for any damage arising from improper use or unauthorised modifications. All risk is borne solely by the device owner. The mixing system must only be started after you have assured yourself that all safety devices are properly installed and fully functional.

Danger

Products which can give off explosive dust/air mixtures must not be processed with this mixing system! The mixing system does not meet ATEX standards! The TOPITEC® AUTOMATIC is not designed for operation in potentially explosive atmospheres.

Proper use of the equipment includes strict compliance with the operating, maintenance and repair instructions of the manufacturer.

The projected service life of the mixing system is approx. 10 years. After 10 years, the device must undergo a complete service by an authorised specialist firm.

2.3 Technical data

2.3.1 Product-specific data

Jar sizes: Jars up to max. diameter 118 mm
Single-patient containers, single-patient jars, jars for stock preparation

2.3.2 Dimensions and weight

TOPITEC® AUTOMATIC (device dimensions) including touch screen

Width: approx. 240 mm
Depth: approx. 240 mm
Height: approx. 680 mm
Total weight: approx. 18 kg

2.3.3 Power supply / electrical parameters

Operating voltage: 1 x 230 V_{AC}
±10%
Phases: 1 Ph / PE
Protective earth conductor: PE (yellow/green)
Frequency: 50 Hz ± 1%
Protection rating: I
Overvoltage category: 2
EMC: Laboratories, domestic environments, SMEs
Protection class: IP 20
Installation standard: DIN EN 61010-1: 2011

Power supply

Fuses: 1x 10A slow-acting and 1x 1A slow-acting
Rated power: approx. 400 W

Warning

The device must be connected to a safety socket. The mixing system is equipped with a fine protection fuse. Medium and general protection must be provided as part of the power distribution system within the building.

The device is protected by 2 fuses (1x 10A slow-acting and 1x 1A slow-acting). The fuse holder is located on the back. At the location of operation, it must be protected by a 10A circuit breaker (characteristic B). When replacing a fuse, use one of the same type, rating and characteristic.

2.3.4 General data

Operating temperature range:

Lower limit temperature:	+ 15 °C
Upper limit temperature:	+ 35 °C
Air humidity:	40%, short-term max. 80%, no condensation
Max. altitude:	≤ 2.000 m above sea level

Storage conditions:

Lower limit temperature:	+ 5° C
Upper limit temperature:	+ 50° C
Air humidity:	40%, short-term max. 80%, no condensation

Danger

Not suitable for operation in potentially explosive atmospheres.

Noise emission:

Sound pressure level:

In accordance with measurement:	$L_{pA} = 55-63 \text{ dB}_{(A)}$
Measurement uncertainty (L_{pA})	2 dB

Warning

Risk from lightning: The power supply to the device must be protected against surges (medium and general protection). Depending on the lightning protection installed on site, it might be necessary to install additional fine protection.

Recommended lighting:

(workplace lighting, to be provided by device owner, according to ASR A3.4, chapter 7.6). Recommendation:
QA workplaces:

$E_x = 500 \text{ lux}$

$E_x = 500 \text{ lux}$

Housing colours:

TOPITEC® AUTOMATIC blue, order no. 025620

TOPITEC® AUTOMATIC white, order no. 025621

3. SAFETY

3.1 Safety instructions and notes

Danger

"DANGER" highlights dangerous situations. Such situations must be avoided! Non-compliance with the instructions might result in serious or even fatal injury.

Warning

"WARNING" highlights potentially dangerous situations. Such situations must be avoided! Non-compliance with the instructions might result in serious or even fatal injury.

Caution

"CAUTION" in conjunction with a warning sign indicates a potentially dangerous situation. Such situations must be avoided! Non-compliance with the instructions might result in minor injury.

Note

"NOTE" highlights instructions that should be adhered to in order to prevent injury. Always follow these instructions to prevent injury, damage to equipment and inconvenience!

Note

Notes in this operating manual / documentation are marked with a "Book" symbol.

3.1.1 Safety symbols in this document

Danger

Danger from electric current

Danger

Danger from running machine (entanglement, crushing)

Danger

Risk of crushing
injury to hands

Warning

When working in areas labelled with this symbol,
wear suitable **protective gloves**.

Warning

When working in areas labelled with this symbol,
wear suitable **protective clothing**.

Warning

Binding instruction: Wear **safety goggles / eye protection**.
Risk of injury to eyes.

Warning

Binding instruction: Wear hairnet.
Observe the hygiene instructions.

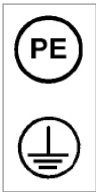
Warning

Risk of **injury from vapours** or dust. During maintenance and repair work,
there is a risk of vapours and dust!

Warning

When working in areas labelled with this symbol,
use suitable **skin protection**.

Note

This symbol indicates the connecting points of the protective earth conductor.

Note

This symbol highlights instructions relating to the protection of the environment.

3.1.2 Device type / identification

Note

The information in this operating manual applies only to devices bearing the type number printed on the cover page. The type plate with the type number is attached to the rear of the mixing system.

When contacting the manufacturer for advice, always quote the following:

- Designation of TOPITEC® AUTOMATIC
- S/N; machine no. 27XXX-01

This ensures that your query can be dealt with quickly.

3.2 Built-in safety systems

The built-in safety devices must be regularly checked and tested by an electrician as described in the table below.

Interval	Method
a = annually	V = visual inspection F = function test M = measurement

Table 1 Test intervals for safety systems

1. Main switch (power switch)

By actuating the main switch, the TOPITEC® AUTOMATIC is connected to / disconnected from the power mains. The main switch is located at the mains filter.

Check	
Interval	Method
a	F

Danger

Before carrying out cleaning, maintenance or repair work, switch off the main switch and also disconnect the power cord from the mains socket.

2. Power supply

The TOPITEC® AUTOMATIC power supply system inside the device consists of a 3 conductor, 1 phase system with a GREEN/YELLOW PE conductor.

Check	
Interval	Method
a	V, M

3. Earthing

The earthing of the electrically conductive components of the TOPITEC® AUTOMATIC must be checked at regular intervals.

Check	
Interval	Method
a	V, M

Warning

The device owner must ensure that unauthorised persons (i.e. persons who are not authorised to operate or service the device) are prevented access to the device.

Operating and maintenance personnel (instructed persons) must be trained and instructed at the location of operation of the TOPITEC® AUTOMATIC. If you have any queries or are unsure how to perform certain tasks, contact your dealer.

Danger

The TOPITEC® AUTOMATIC contains internal components that are electrically powered. Its internal parts do not need to be serviced and there is no need to open the housing of the device! Please ensure that the TOPITEC® AUTOMATIC is only opened by an electrician for maintenance purposes and otherwise remains closed!

Warning

This operating manual forms an integral part of the scope of delivery of the mixing system. A copy of this manual must be accessible at all times to operating and maintenance personnel.

All safety instructions in this document must be strictly adhered to.

When selling the device, this manual must be handed over to the new owner.

Warning

Never disable or otherwise interfere with the safety devices of the mixing system. After a failure, repair or service, you must ensure that all safety devices are correctly installed and working properly.

3.3 Safety precautions (to be implemented by device owner)

Please note that the device owner is responsible for the proper instruction of all operating and maintenance personnel as regards:

- the **safety devices** of the TOPITEC® AUTOMATIC;
- compliance with all safety instructions in this manual.

Compliance must be monitored. A copy of this operating manual must be filed safely for future reference. All inspection and maintenance intervals must be strictly observed.

The various sections of this operating manual contain instructions for different groups of personnel:

- The chapters referring to operation and operating modes are **written for** instructed personnel.

The instructions in the chapters referring to **troubleshooting and repair** are **aimed at qualified electricians**, as all tasks described in these chapters must only be performed by qualified electricians.

Instructed person (operator)

A person who has been instructed by a qualified electrician in the proper operation of the device, the associated risks and preventive measures, the purpose and functions of the safety devices and best safety practice.

Electrician

A person (maintenance/repair technician) who, due to his/her professional training and experience is in a position to identify and prevent risks.

3.4 Duties of device owner

Note

In EFTA countries, the relevant national regulations based on framework directive 89/391/EEC and the associated specific directives, in particular 2009/104/EC regarding the minimum health and safety precautions in connection with the operation of equipment by employees, as amended, apply and must be strictly adhered to.

In Germany, the Ordinance on Industrial Safety and Health of 03/02/2015 (BetrSichV) applies. Also observe the German Ordinance on the Operation of Pharmacies ApBetrO and the BAK Guidelines.

The device owner is responsible for compliance with all relevant statutory regulations in connection with

- occupational health and safety (accident prevention regulations, etc.), for example by devising internal operating instructions, providing personal protective equipment (PPE) and arranging for preventive check-ups;
- the safety of the equipment (e.g. protective equipment, work instructions, process risk assessments, and maintenance schedules);
- the handling of hazardous substances;
- the procurement of materials (material safety data sheets, hazardous substances lists);
- product disposal (waste disposal regulations);
- the disposal of the device (decommissioning, waste disposal regulations);
- cleaning (cleaning agents and their disposal);
- the protection of the environment;
- compliance with the ApBetrO, the guidelines published by the German Federal Chamber of Pharmacists BAK, known as the BAK Guidelines, and all other relevant statutory regulations.

The device owner is also obliged to

- provide and regularly check the personal protective equipment (PPE): lab coats, protective gloves, safety goggles, hoods/hairnets, etc.
- carry out regular risk assessments of the individual workplaces, including observation of hazardous substances and fire risks;
- issue binding operating instructions and train personnel in the safe operation of the device (training of operating personnel).
- issue work instructions for staff.
- enforce strict rules to ensure that personnel do not operate the device under the influence of alcohol or other drugs, or when extremely tired.
- implement procedures that ensure that all safety-relevant faults of the device or its components are reported without delay to the supervisor.
- devise a skin protection and hygiene plan.

The device owner is also obliged to

- provide suitable safety equipment such as hand-held fire extinguishers in the required quantity and of suitable size and type at easily accessible locations near the device;
- protect the TOPITEC® AUTOMATIC through a potential equalisation system that conforms to the statutory requirements (in Germany: Ordinance on Industrial Safety and Health);
- provide adequate lighting at the workplaces (conforming to ASR A3.4);
- ensure that the ventilation system at the workplaces remains switched on outside working hours to maintain a minimum air flow at all times (elimination of germs);
- regularly inspect the working area to ensure that tripping hazards are eliminated or suitably marked.

Connections:

Prior to operating the TOPITEC® AUTOMATIC it is essential that the device owner ensures that the connected loads and the local regulations (e.g. for the electrical connection) are observed.

Note

Lighting:

The device owner must provide adequate and uniform lighting of the prescribed illumination rating in the area of the TOPITEC® AUTOMATIC. Recommended: **500 lux** (QA workplaces **500 lux**), depending on site conditions (maintenance level; in Germany: ASR A3.4).

Note

Noise:

The TOPITEC® AUTOMATIC does not produce noise at a damaging sound pressure level (for limits, see German Noise and Vibration Protection Ordinance LärmVibrArbSchV). **Sound level > 85 / 80 dB_(A)**

Warning

At least one **safety training session** per year (and also upon appointment of a new employee), covering all risks and safety measures, must be provided for all staff members, based on this operating manual and in a language that is easily understood by staff. Completion of the training must be confirmed by signature.

3.5 Safety tests

performed at the factory by the device manufacturer

1. Risk analysis based on Machinery Directive 2006/42/EC (Annex I) and DIN EN ISO 12100:2011.
2. Sound pressure measurement (no report required, as LPA < 70 dB(A)) according to Machinery Directive, Annex I
3. Inspection and testing according to DIN EN 61010-1: 2011 2011

3.6 Safety checks

The safety tests and checks of the TOPITEC® AUTOMATIC have been performed by the manufacturer at the factory. All relevant procedures and results have been documented. The device owner is obliged to carry out repeat safety checks and tests according to DGUV A3.

4. GENERAL SAFETY INSTRUCTIONS

4.1 Dangers

Observe the information regarding the safety systems, and the safety notes in this operating manual. The device is controlled through the keypad. During operation, keep the area around the TOPITEC® AUTOMATIC free of clutter to ensure unobstructed access.

Danger

When carrying out maintenance or repair work, there is a risk of injury from crushing!

Danger

Pay attention to all moving parts while working. Risk of entanglement! When working in the danger area, wear tight-fitting clothing (no necklaces, rings, ties, long hair, long earrings, etc.). Do not wear protective gloves (risk of entanglement).

Danger

When carrying out maintenance or repair work, avoid contact with **electrically powered parts!**

Protect the housing against the ingress of water or other foreign substances. Do not place objects on the device, as this could lead to overheating.

4.2 Danger areas of TOPITEC® AUTOMATIC

For operating personnel, the operating area includes the keypad and the other control elements as well as the areas that need to be accessed to set up the TOPITEC®AUTOMATIC and to clean it. When carrying out maintenance or repair work, the entire area around the TOPITEC® AUTOMATIC **becomes a danger area** that must only be accessed by electricians who must strictly adhere to the safety instructions.

4.3 Operating and maintenance personnel

Operating and maintenance personnel include all persons involved in the transportation, installation, operation and cleaning of the TOPITEC® AUTOMATIC. (For responsibilities and authorisations, see page 18).

1. The TOPITEC® AUTOMATIC must only be operated by suitably trained and authorised personnel.
2. The device owner must ensure that there are clear instructions regarding the operation of the TOPITEC® AUTOMATIC in place, and that all personnel are aware of their responsibilities so as to avoid confusion that might result in a dangerous situation.
3. For all work with and on the device (operation, cleaning, maintenance, repair, etc.), personnel must strictly adhere to the shut-down procedures described in this operating manual.
4. Operating personnel must refrain from any action that might compromise the safety of the TOPITEC® AUTOMATIC.
5. The device owner must take suitable measures to ensure that only authorised personnel have access to the TOPITEC® AUTOMATIC.
6. Operating personnel are obliged to notify their supervisor without delay, if they observe any changes to the TOPITEC® AUTOMATIC that might compromise its safety.
7. The device owner must ensure that the TOPITEC® AUTOMATIC is always in proper working order.
8. Operating personnel must comply with all statutory regulations and observe the specific instructions for the substances to be processed. Information regarding proper hygiene procedures can be found in the ApBetrO, the hygiene guidelines published by Gesellschaft für Dermopharmazie GD for the preparation of sterile pharmaceutical products, the BAK Guidelines, statutory regulations, etc.
9. The device owner must ensure that all personnel wear the required personal protective equipment. Compliance must be monitored.

Danger

PPE:

Wear tight-fitting clothing, preferably without sleeves. Do not wear scarfs, ties, necklaces, rings, etc. Cover your hair with a hairnet. Do not operate the device with gloves, as there is a risk of entanglement at the shaft.

Wear safety goggles to protect your eyes against harmful substances from ointments, creams, powders, etc. While the machine is in operation, do not reach into its mechanical parts. Do not touch the shaft, as this could lead to serious injury from crushing or shearing.

Moving the safety bracket (see also page 32) upwards leads to the device shutting down. After an emergency-stop, the device must be restarted.

Information regarding proper hygiene procedures can be found in the Ap-BetrO, the hygiene guidelines published by Gesellschaft für Dermopharmazie GD for the preparation of non-sterile pharmaceutical products (known as GD Hygiene Guidelines), the BAK Guidelines, statutory regulations, etc.

4.4 Replacement of parts

Please note that third-party replacement or wear parts have not been tested or approved by us. The installation and/or use of such products might therefore impair the operation of the TOPITEC® AUTOMATIC. We shall not accept any liability for damage caused by the use of third-party parts or accessories.

Warning

After the replacement of wear parts, check all safety devices for proper functioning.

Do not start the mixing system before you have assured yourself that all safety devices are working properly.

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below:

Danger

Prior to any cleaning, maintenance or repair tasks (to be performed by an electrician), shut down the device as described

1. Lower the slide to its initial position and remove the jar.

2. Disconnect the TOPITEC® AUTOMATIC from the power supply:

- Set the main switch (power switch) to position "0".
- Secure the TOPITEC® AUTOMATIC against inadvertent restart by disconnecting the power cable from the socket.

2.1 For maintenance and repair work only:

- Check the device to ensure that it is fully de-energized.
- Cover or otherwise insulate adjacent parts that remain powered.

Danger

Non-compliance can result in serious or even fatal injury (from accidental voltage transfer)!

Capacitors might retain some dangerous residual voltage.

Before commencing work, check the capacitors and de-energize them slowly.

If a task cannot be performed when the device is shut down as described above, contact your dealer or carry out the task with due regard to the internal risk assessment.

5. STANDARD EQUIPMENT / INSTALLATION / COMMISSIONING

5.1 Scope of delivery of TOPITEC® AUTOMATIC

The delivery includes:

1. TOPITEC® AUTOMATIC
2. Operating manual
3. Standard equipment

5.2 Basic equipment

THE TOPITEC® AUTOMATIC

is available in two versions

Order no. 025620 - Housing colour blue

Order no. 025621 - Housing colour white

Power cable, order no. 080850

TOPITEC® holder for jar 20/30 g

for TOPITEC® jar sizes 20 g and 30 g, order no. 026631

TOPITEC® holder for jar 50-200 g

for TOPITEC® jar sizes 50 g to 200 g, order no. 026632

Standard kit of TOPITEC® single patient jars, order no. 025688

- 1 mixing tool for large containers
- 2 mixing lids for large containers
- 6 sealing rings for mixing lid

Tool shaft for mixing discs,

TiN-plated, order no. 025680

Receipt booklet

including TOPITEC® jars user instructions, order no. 038053

Fig. 1-2 Colours of TOPITEC® AUTOMATIC

5.3 Installation

5.3.1 Transport and packaging

All mixing systems are carefully inspected before they are packed. However, the equipment might become damaged during transport.

5.3.1.1 Delivery (also for spare and replacement parts)

Incoming inspection

- Upon receipt of the delivery, check the goods for completeness, referring to the delivery docket!

In the event of damage:

- Check the transport packaging for visible damage!
- Check the goods for visible damage.

TOPITEC® AUTOMATIC packaging

1. Place the cardboard box on the floor, observing the "this side up" mark.
2. Cut the top of the packaging and fold down the four cover flaps.
3. Remove all packaging as far as this is possible at this stage.
4. Use two people to lift the device (weight approx. 18 kg) out of the box and place it at the designated spot.
5. Remove all remaining packaging.

Complaints procedure

If the goods have been damaged during transport:

- Immediately contact your dealer!
- Keep the packaging (for further inspection by the shipping company or for return of the goods).

Packaging for return:

Use the original packaging.

If the original packaging is not available:

- Contact our customer service and request a packaging kit.
- If you have any queries regarding the packaging and securing of the goods for transport, contact your dealer.

5.3.1.2 Temporary storage

The transport packaging of the TOPITEC® AUTOMATIC and spare parts is designed for storage for maximum 3 months.

Note

Store the TOPITEC® AUTOMATIC in a dry room.

Storage conditions:

- See chapter 2.3.4 "Technical data".

Warning

Never clean the TOPITEC® AUTOMATIC with a steam cleaner or a water jet. Water and dirt might otherwise enter the mixing system, causing damage.

5.4 Transport

5.4.1 Technical transport instructions

Warning

When transporting the device, the **weight** and the **centre of gravity** of the TOPITEC® AUTOMATIC must be observed (see Table 2). During transport, the TOPITEC® AUTOMATIC may tip over. To prevent this, always observe its **centre of gravity**. Secure the packaging of the TOPITEC® AUTOMATIC with appropriate slings **prior to transport if necessary**.

When **transporting the TOPITEC® AUTOMATIC**, the possible tipping over of system parts must also be taken into consideration when determining the space requirements (height of the components).

Assembly / transport unit	Weight	Centre of gravity	Lashing points	Lifting gear
TOPITEC®AUTOMATIC (heaviest transport unit)	approx. 18 kg (net)	Top-heavy	./.	by hand (1 person)

Table 2 Weights of transport units

5.5 Installation and assembly

Warning

Install the device on a firm and stable, non-slipping surface. The base of the device must be fully supported and positioned at least 20 cm away from the front and rear edges of the worktop, and 50 cm from the right and left edges of the worktop. Above the device, there must be a clearance of at least 100 cm to ensure proper ventilation and fire safety. Do not position the device in a highly frequented area or corridor. The device owner must ensure that the workplace of the device meets the general health & safety standards.

5.6 Commissioning

5.6.1 Start-up

Remove all packaging and protective material before setting up the TOPITEC® AUTOMATIC (see chapter 5.3.1.1). The TOPITEC® AUTOMATIC must be placed on a level and firm surface. When choosing the location of installation, observe the information in section 5.5 (Installation and assembly).

The TOPITEC® AUTOMATIC is not designed for operation in potentially explosive atmospheres. When handling and using hazardous substances, always strictly adhere to the relevant safety requirements. For more information, refer to the BAK Guidelines and the German Hazardous Substances Ordinance.

- Connect the power cord to the mains filter socket at the rear of the device.
- Connect the plug of the power cable to the mains socket. Ensure that the power socket is correctly earthed!
- Switch on the device, following the instructions in chapter 6.

The power cable socket of the TOPITEC® AUTOMATIC is located at the rear of the device. Compare the mains voltage with the rated voltage of the system (230 V). The mains plug should always be easily accessible.

6. OPERATION

Warning

The TOPITEC® AUTOMATIC must only be operated by instructed personnel who have been trained in its functions and control.

Note

Before starting production, ensure that **there are no objects** (e.g. tools, residual material, etc.) on or near the TOPITEC® AUTOMATIC mixing system.

6.1 Connection to power supply, interfaces

- 1. RS 232 interface
- 2. Fuse holder for slow-acting 1A fuse
- 3. Power switch (main switch)
- 4. Fuse holder for slow-acting 10A fuse
- 5. Power socket (3-pin)

Fig. 1-3 Power connection, interfaces, fuse holders

The rear panel of the TOPITEC® AUTOMATIC features the following control devices and ports:

- 1. **RS 232 interface**
for the printer connection (see 6.2)
- 2. **Fuse holder (recessed)** containing 1 slow-acting 1A fuse.
- 3. **Main switch** (power switch)

When this switch is ON, the TOPITEC® AUTOMATIC is powered from the power mains.

Warning

If the main switch needs to be switched off for cleaning, maintenance or repair work, always disconnect the power cable from the power socket (to prevent inadvertent switching on).

- 4. **Fuse holder (recessed)** containing 1 slow-acting 10A fuse.
- 5. **Power socket (3-pin)**
with phases including PE conductor; socket for IEC-60320 C13/C14 cable plug. The power cable must be connected to a properly earthed power socket with a PE conductor.

6.2 Interfaces of TOPITEC® AUTOMATIC

For the documentation of your preparations and production runs, we recommend connecting a SARTORIUS data printer (model YDP20-0CE) to your TOPITEC® AUTOMATIC

1.

1. First check whether the printer is correctly configured for connection and printing. The baud rate must be set to 9600. For detailed instructions, please refer to the user manual of the printer.

2. Connecting the printer

Connect the printer power cable to the socket. The printer is automatically switched on. The display shows the time and the activated numerator (NUM). Deactivate the numerator by pressing the [N/N] button.

3.

3. The serial port for PC connection (RS 232) is located at the rear of the TOPITEC® AUTOMATIC.

Connect the serial cable (prod. no. 25 025663) to the serial port at the rear of the printer (labelled DATA).

4.

4. The printer cable can be connected.

Ensure that all adapters are properly inserted in the ports. The printer can be connected before the TOPITEC® AUTOMATIC is started or during production.

6.3 Safety bracket

When the safety bracket is touched, the mixing system is instantly stopped.

The safety bracket prevents injury from being caught between the tool shaft adapter and the jar during production. If the safety bracket is pushed up while the slide is moving upwards, the mixing process is immediately stopped.

Safety bracket

Fig. 1-4 Safety bracket

6.4 Control elements of TOPITEC® AUTOMATIC

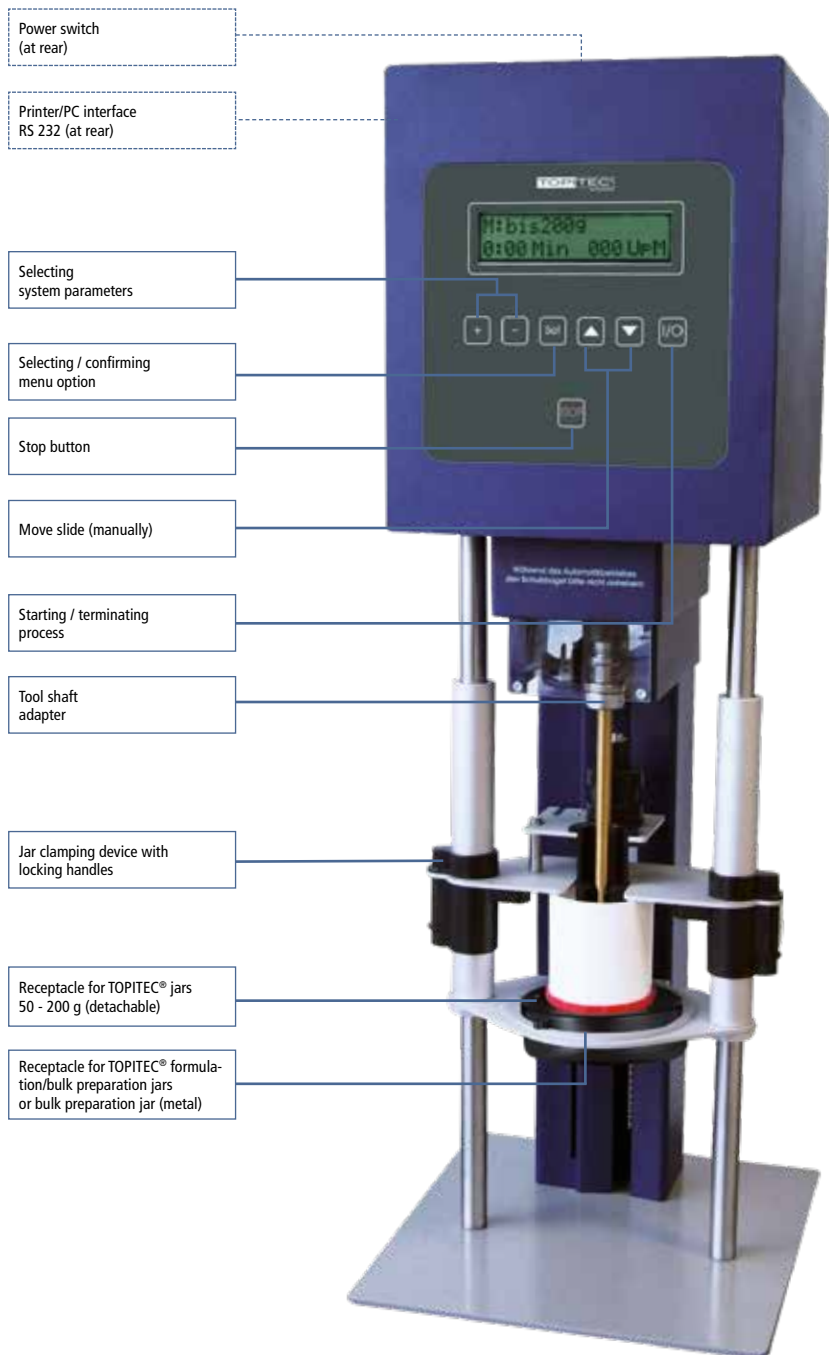
Fig. 1-1 Control elements of TOPITEC® AUTOMATIC

The control of the TOPITEC® AUTOMATIC mixing system occurs via a keypad in the form of symbols and button functions. Do not apply force when touching the surface. If the keypad buttons are touched with an implement, the keypad might become dirty or damaged.

6.5 Operating modes and control

6.5.1 Manual mode for cleaning

Activate the manual slide function to clean the TOPITEC® AUTOMATIC. You can now move the slides in any direction you wish with the arrows.

Note

Before starting production, ensure that there are **no** objects (e.g. tools, residual material, etc.) on or near the TOPITEC® AUTOMATIC.

6.5.2 Automatic mode

Automatic mode is the standard mode for production. Enter the formulation and start the process. Once the stirring process is complete and following subsequent manual unlocking of the tool shaft, it is possible to move the slide back to its home position via the "Arrow down" button.

6.5.3 Switching on the mixing system

Connect the TOPITEC® AUTOMATIC with the supplied power cable to a properly earthed power socket. Connect the power switch.

6.5.4 Operation

Warning

- Before starting the device, check the power cable and the main switch for proper connection/functioning.
- Switch on the main switch. The device starts up.
- The TOPITEC® AUTOMATIC mixing system is controlled through the keypad.

Warning

When in operation, the device must be constantly supervised.

6.5.5 Switching off the mixing system

- After the mixing process is completed, remove the mixing jar from the holder.
- Switch off the main switch.
- Clean the TOPITEC® AUTOMATIC and its accessories.
- Switch off the power switch and disconnect the power plug from the socket if the device is not required.

6.5.6 Restarting mixing system after error or fault

If the mixing system has been affected by a fault or error, first have the problem eliminated by a specialist technician. After a thorough inspection of the device, restart it.

6.5.7 Restarting after power failure

When the power returns, move the mixing system by hand to its initial position and check it. If everything is OK, set the device to automatic mode.

MODE OPTIONS / BUTTON FUNCTIONS

- **M:** to 200 g
- **M:** from 300 g
- **Comminution**
- **Powder mixing**

Mode selection for once-off production of single preparation

Press 1 x **Set**: "M: to 200 g" flashes

Select with **+/-**: "Mode"

Press 1 x **Set**: "Min" flashes

Set with **+/-**: "Min"

Press 1 x **Set**: "rpm" flashes

Set **+/-**: "rpm" with

Press **Set** 1x to confirm entry

Press 1 x **I/O** to start the process

Slide moves up & tool shaft engages.
Programme is started ("rpm" flashes; if necessary, adjust rpm by pressing **+/-**).
Program time countdown begins.

"Program terminated"
"Release tool shaft"
"Move slide"

Lift the safety bracket and keep it upwards, unlock tool shaft from the shaft retainer, pressing the **▼** button until the slide has moved downwards with the jar clamped in place.

Automatic: Default settings

Button functions

- +/-** Set values for "min.", "rpm"; select mode, or "Prep. no."
- Press 1x "Manual slide movement" to move slide based on default settings (M: to 200 g, 0:00 min., 000 rpm)
- Set** Select menu option (mode/min./rpm/pause/prep. no.) and confirm
- ▲** Actuate when prompted by system: "Release tool shaft" and "Move slide". Do not forget to release the tool shaft!
Lift the safety bracket!
- I/O** When the container is engaged, the next cycle is started.
Alternatively, the process is terminated prematurely. Next the stirring programme is continued following a safety bracket interruption.
- Stop** Only press this button if there is an error or malfunction! (After switching off the TOPITEC® AUTOMATIC system it must be restarted, see TOPITEC® manual, section "Troubleshooting") Exception: Change from save mode to default mode.

TOPITEC® AUTOMATIC QUICK START

Initial position:
Mixing container clamped in slide Default settings:
M: to 200 g, 0:00 min. , 000 rpm

Setup for once-off preparation M: to 200g

Press 1 x **Set**: "M: to 200 g" flashes

Press 1 x **Set**: "Min" flashes

Set with **+/-** "Min"

Press 1 x **Set**: "rpm" flashes

Set "rpm" with **+/-**

Press **Set** 1x to confirm entry

Press 1 x **I/O** to start the process

Slide moves up & tool shaft engages.

Programme starts ("rpm" flashes and could be changed if necessary by pressing **+/-**).

Program time countdown begins.

"Program terminated"
"Release tool shaft"
"Move slide"

Lift safety bracket and keep it in the upwards position, release tool shaft from shaft retainer, pressing the **▼** button down until the slide has moved downwards with the jar clamped in place.

Automatic: Default settings

Single-step preparation

Press **Set** 1x: "Prep. no." (e.g. 1/1) flashes

Press **I/O** 1x to start process

Slide moves up & tool shaft engages.

Program time countdown begins.

"Program terminated"
"Release tool shaft"
"Move slide"

Lift safety bracket and keep it in the upwards position, release tool shaft from shaft retainer, pressing the **▼** button down until the slide has moved downwards with the jar clamped in place.

Press **Stop** 1x: Default settings

Button functions

+/- Set values for "min.", "rpm"; select mode, or "Prep. no."

- Press 1x "Manual slide movement" to move slide based on default settings (M: to 200 g, 0:00 min., 000 rpm)

Set Select menu option (mode/min./rpm/pause/prep. no.) and confirm

▼ Actuate when prompted by system: "Release tool shaft" and "Move slide". Do not forget to release the tool shaft! Lift the safety bracket!

I/O When the container is engaged, the next cycle is started. Alternatively, the process is terminated prematurely. Next the stirring programme is continued following a safety bracket interruption.

Stop Only press this button if there is an error or malfunction! (After switching off the TOPITEC® AUTOMATIC system it must be restarted, see TOPITEC® manual, section "Troubleshooting"). Exception: Change from save mode to default mode.

Define single-step or two-step preparations (max. 3 programming steps) M: to 200g

Press **Set** 1x: "Prep. no." flashes

Select "Prep. no." (e.g. 1) with **+/-**

Press **Set** 1x: "M: to 200 g" flashes

Press **Set** 1x: "Min" flashes

Set "min." with **+/-**

Press **Set** 1x: "rpm" flashes

Set "rpm" with **+/-**

Press **Set** 1x: "pause" flashes

Set "pause" with **+/-**: "Min" flashes OR
press **Set** to skip "pause": "min." flashes

Two-step preparations

Set "min." with **+/-**

Press **Set** 1x: "rpm" flashes

Set "rpm" with **+/-**

Press **Set** 1x: "pause" flashes

Set "pause" with **+/-**: "Min" flashes OR
press **Set** to skip "pause": "Min" flashes

Press **Set** 1x: "Prep. no." (e.g. 1/2) flashes

Press **I/O** 1x to start process

Slide moves up & tool shaft engages.

Program time countdown begins.

"Program terminated"
"Release tool shaft"
"Move slide"

Lift safety bracket and keep it in the upwards position, release tool shaft from shaft retainer, pressing the **(V)** button down until the slide has moved downwards with the jar clamped in place.

Press **Stop** 1x: Default settings

Three-step preparations

Set "min." with **+/-**

Press **Set** 1x: "rpm" flashes

Set "rpm" with **+/-**

Press **Set** 1x: "pause" flashes

Set "pause" with **+/-**: "Min" flashes OR
press **Set** to skip "pause": "Min" flashes

Set "min." with **+/-**

Press **Set** 1x: "rpm" flashes

Set "rpm" with **+/-**

Press **Set** 1x: "Prep. no." (e.g. 1/3) flashes

Press **I/O** 1x to start process

Slide moves up & tool shaft engages.

Program time countdown begins.

"Program terminated"
"Release tool shaft"
"Move slide"

Lift safety bracket and keep it in the upwards position, release tool shaft from shaft retainer, pressing the **(V)** button down until the slide has moved downwards with the jar clamped in place.

Press **Stop** 1x: Default settings

MIXING MODE / PREPARATION SETUP (ONCE-OFF)

1.

1. If you wish to process a formulation that you do not require at any later stage, you have the option to adjust the default settings according to your requirements without the need to save these settings. To do this, press **(Set)** 2x. "min." flashes on the display, indicating that you can now adjust the time. To do this, press the **(+/-)** buttons as required. Press the **(Set)** button. "rpm" flashes on the display, indicating that you can now adjust the speed.

To start production, press the **(I/O)** button.

2.

2. The slide moves to its end position, and the message to the left is shown on the display:

The TOPITEC® AUTOMATIC guides the tool shaft into the drive shaft sleeve automatically and locks it in position. After a short run-up time of the torque motor, production starts with the user-defined settings.

The speed and time remaining until the end of the mixing process is shown on the display.

The TOPITEC® AUTOMATIC display also shows a number (1 to 99) referred to as the production number. It is generated automatically and shown during the entire production time in the top right corner of the display (here: 11).

3.

3. "rpm" flashes on the display, indicating that you can adjust the speed throughout the mixing process by means of the **(+)** / **(-)** buttons.

MIXING MODE / PREPARATION SETUP (ONCE-OFF)

4a.

- 4a. After completion of the mixing process, message "Process terminated" is displayed.

4b.

- 4b. After approx. 10 seconds, an acoustic signal is generated, and the display changes to "Release tool shaft".

5.

5. Using one hand, lift the safety bracket to the stop, pulling the lock ring of the shaft sleeve upwards. With the second hand, press the  button. The slide returns to its initial position. Release the two locking handles at the sides and pull the locking device upwards. The TOPITEC® jar can now be removed from the device!

WORKING WITH CUSTOMISED SETTINGS

The machine's settings for mixing can of course be modified by adjusting the values for "min." and "rpm". In the example below, the following parameter values are entered:
M: to 200 g; 1:30 min.;1500 rpm.

1. Display with default settings
Press **Set** 2x.

1.

2. Press **+** to set the time to 1:30 minutes

2.

3. Press **Set** 2x.
Press **+** to change the speed (rpm)

3.

4. To start the process, press **I/O** button.

4.

SEQUENCE OF STORED PREPARATIONS

Switch on the system... > M: to 200 g
0:00 min. 000 rpm

Press **+** to select... > Prep. no. 1 to prep. no. 99

Press **-** to select... > move slide

Setting up single-step preparations

If you wish to produce the same preparation more than once, you can save the respective system settings in one of the 99 available memory slots so that they can later be called up at the push of a button. After switching on the system, the default settings are displayed.

MIXING MODE / REPRODUCIBLE PREPARATION SET-UP

Setting up single-step preparations

1.

1. If you wish to produce the same preparation more than once, you can save the respective system settings in one of the 99 available memory slots so that they can later be called up at the push of a button. After switching on the system, the default settings are displayed.

2.

2. To call up a memory slot, press **+** (here: 1x).
"prep. no." flashes on the display.

3.

3. Press **Set** :
"M: to 200g" (mode display) flashes. Press **Set** again:
"M: to 200g" is continuously shown. "min." flashes.

4.

4. Using the **+** button, enter the desired time in minutes and seconds. If you have inadvertently selected a value that is too high, press the **-** button to correct it. Press **Set** 1x.
You can now adjust the speed.

5.

5. Using the **+** button, enter the desired speed.

6.

6. Now press the **Set** button 3x. The preparation number is shown on the display:

7. All entered values are saved under the respective prep. no. (here: 1). In future you only need to call up the "Prep. no. 1" and production can start. To start production, press the **I/O** button. Alternatively, you can set other preparations in the system or press **Stop** to return to the TOPITEC® AUTOMATIC default settings.

MIXING MODE / REPRODUCIBLE PREPARATION SET-UP

Setting up multiple-step preparations

Each preparation can be divided into three separate steps with individual time (min.), speed (rpm) and pause settings.

Example:

- "Prep. no. 10"
- step 1: 0:30 minutes at 500 rpm, pause time 1:00 minute
 - step 2: 2:00 minutes at 1500 rpm, pause time 2:30 minutes
 - step 3: 1:30 minutes at 2500 rpm

1.

1. Press **+** to select "Prep. no. 10"

2.

2. Press **Set** 2x:
"min." flashes.

3.

3. Using the **+** button to enter the time in minutes (here: 0:30). Press the **Set** button again:
"rpm" flashes. Using the **+** button to enter the speed (here: 500). Confirm with **Set**.

"pause" flashes.

4.

4. Using the **+** button, enter the pause time (here: 1:00 and confirm with **Set**).

"min." flashes.

MIXING MODE / REPRODUCIBLE PREPARATION SET-UP

5.

5. You are now ready to enter the second step of preparation. Repeat the above steps (here, 2:00 min.; 1500 rpm; 2:30 pause

"pause" flashes.

6.

6. Confirm with the **Set** button. You are now ready to enter the settings for step 3 of the preparation.

"min." flashes.

7. Using the **+** and **Set** button, enter the relevant values for time and speed (here: 1:30 min.; 2500 rpm). Confirm the rpm entry with **Set**. In our example the preparation is saved in the TOPITEC® AUTOMATIC system under "prep. no. 10".

You can now start production by pressing **I/O** (TOPITEC® jar in place?), set other preparations in the system or press **Stop** to return to the TOPITEC® AUTOMATIC default settings.

Preparations can consist of one, two or three steps. To configure a preparation with two steps, follow the instructions above and confirm your entries by pressing **Set** twice after entry of the second "rpm" value.

PAUSE FUNCTION

1.

1. For multiple-step preparations, you have the option to define pause times (max. 30 minutes per step).

This enables you to include a cooling phase in the process.

The pause time is automatically added by the TOPITEC® AUTOMATIC to the overall production time and shown on the display during production.

To prevent inhomogeneous results, the TOPITEC® AUTOMATIC stirs the product 3 times every 2 minutes at low speed (300 rpm) during the pause.

CHANGING VALUES IN SAVED PREPARATIONS

1.

1. Press **+** to call up the respective preparation.
(Example: Prep. no. 10; 1. step NEW: 1:20 minutes at 2000 rpm).

"Prep. no." flashes.

2.

2. Press **Set** 2x. The time entered for the 1st step is displayed.
Press **+** to change the time to 1:20 minutes. Press **Set** again.

"rpm" flashes.

Press **+** to change the value to 2000 rpm.

Repeatedly press **Set** until "prep. no." is displayed again (flashing).

The preparation parameters are now changed and saved in the TOPITEC® AUTOMATIC.

Press **Stop** to return to the TOPITEC® AUTOMATIC default settings.

DELETING SAVED PREPARATIONS

1.

1. Press **+** to call up the respective preparation.
(Example: Prep. no. 10):

"Prep. no." flashes.

2.

2. Press **Set** 2x. The time entered for the 1st step is displayed
Press the **-** button to set the time to 0:00 and then
press **Set** again.

"Prep. no." flashes.

All additional steps and the set "Prep. no. 10" are deleted or e.g. a memory slot can now be used for a new preparation. Press **Stop** to return to the TOPITEC® AUTOMATIC default settings.

7. PRICING

7.1 aponorm® packaging units for TOPITEC® mixing systems

The latest edition of the "Hilfstage für Apotheken" that applies in Germany contains the agreed prices for the aponorm® jars with metering bottom (TOPITEC® jars). For your convenience, the details of the packing units (aponorm® jars with metering bottom (TOPITEC® jars)), aponorm® tamper-evident jars (TOPITEC® tamper-evident jars), aponorm® single-patient jars (TOPITEC® single-patient jars) and other TOPITEC® accessories are compiled in the table below:

aponorm® jars with metering bottom (TOPITEC® jars)				
Order no.	Colour/design	Capacity	PZN (1 unit)	SU (quant.)
035250	red/white	20 g / 28 ml	0775126	20
035251	red/white	30 g / 42 ml	0775161	20
035252	red/white	50 g / 70 ml	0775190	20
035253	red/white	100 g / 140 ml	0775221	15
035255	red/white	150 g / 200 ml	2381704	10
035254	red/white	200 g / 250 ml	0775244	10
035230	blue/blue	20 g / 28 ml	2684385	20
035231	blue/blue	30 g / 42 ml	2684391	20
035232	blue/blue	50 g / 70 ml	2684416	20
035233	blue/blue	100 g / 140 ml	2684422	15
035235	blue/blue	150 g / 200 ml	0764683	10
035234	blue/blue	200 g / 250 ml	2684439	10
035425	orange/cosmetics	20 g / 28 ml	2726379	20
035426	orange/cosmetics	30 g / 42 ml	2726385	20
035427	orange/cosmetics	50 g / 70 ml	2726391	20
035428	orange/cosmetics	100 g / 140 ml	2726422	15
035429	orange/cosmetics	150 g / 200 ml	2726439	10
035430	orange/cosmetics	200 g / 250 ml	2726445	10

aponorm® tamper-evident jars (TOPITEC® tamper-evident jars) & accessories				
Order no.	Colour/design	Capacity	PZN (1 unit)	SU (quant.)
033150	red/white	20 g / 28 ml	7317326	20
033151	red/white	30 g / 42 ml	7317332	20
033152	red/white	50 g / 70 ml	7317349	20
033153	red/white	100 g / 140 ml	7317355	15
033155	red/white	200 g / 250 ml	7317361	12
096090	Ointment spatula	100 mm	—	20
096092	Ointment spatula	100 mm	—	15
096091	Ointment spatula	150 mm	—	12

INDIVIDUELL

EINZIGARTIG

For jars and ointment spatulas with your own Pharmacy imprint – contact us.

Fig. 1-5 aponorm® packaging units for TOPITEC® mixing systems

aponorm® single-patient jars (TOPITEC® single-patient jars)

Order no.	Colour/design	Capacity	PZN (1 unit)	SU (quant.)
035260	red/white	300 g / 408 ml	0775379	10
035261	red/white	500 g / 690 ml	0775416	10

aponorm® jars for stock preparation (TOPITEC® jars for stock preparation)

Order no.	Colour/design	Capacity	PZN (1 unit)	SU (quant.)
035263	red/white	1,000 g / 1,260 ml	3451051	4

ACCESSORIES

The table below contains information you might require for the pricing of TOPITEC® accessories or for your stock management:

TOPITEC® accessories

Order no.	Item	Content/length	PZN (1 unit)	SU (quant.)
025690	Mixing discs	20 g and 30 g	0775304	50
025691	Mixing discs	50 g to 200 g	0775310	50
035270	Ointment cannula	1.5 mm	0775327	30
035271	Reducer element	2.5 mm	0775333	30
035275	Dosing sleeve	–	0775362	200

These accessories and their prices can be entered into your formulation pricing program. For details, please contact your software supplier.

Metering tips for aponorm® jars with metering bottom (TOPITEC® jars)

made in PP, red, SU 200

order no. 035275; PZN 0775362

Ointment cannula for aponorm® jars with metering bottom (TOPITEC® jars)

for targeted application, e.g. in body cavities,
1.5 mm upper Ø, made in PP, white, SU 30

order no. 035270; PZN 0775327

Ointment cannula for aponorm® jars with metering bottom (TOPITEC® jars)

for accurate dosing of low-viscous formulations,
top Ø 2.5 mm, made in PP, white, SU 30

order no. 035271; PZN 0775333

8. ACCESSORIES FOR TOPITEC® AUTOMATIC

8.1 aponorm® jars (TOPITEC® jars)

Receipt book, incl. TOPITEC® jars
Use instructions
order no. 038053

aponorm® single-patient jars (TOPITEC® single-patient jar
made in polypropylene (a product of the aponorm® packaging range),
with batch certificate, SU 10
300 g / 408 ml single-patient jar, **order no. 035260; PZN 0775379**
500 g / 690 ml single-patient jar, **order no. 035261; PZN 0775416**

Mixing tool for aponorm® single-patient jars
(TOPITEC® single-patient jars)
SU 1, **order no. 025695**

The mixing tool is made in stainless steel (exclusively for the single-patient jar). Its shaft is protected by an extra resistant TIN alloy coat. The mixing element ($\varnothing \times H = 103.5 \times 4.0$ mm) is equipped with a wear-proof plastic (PTFE) coat.

Mixing lid for aponorm® single-patient jars
(TOPITEC® single-patient jar)

The mixing lid is made in white POM. It features a centre guide hole or the tool and is equipped with a recessed silicone O-ring seal.
SU = 2, **order no. 025682 (white)**

Fig. 1-6 Accessories TOPITEC® mixing systems

**aponorm® jar for stock preparation
(TOPITEC® jar for stock preparation)**

for the production and storage of semi-solid formulations produced with TOPITEC® AUTOMATIC II, TOUCH and EXPERT mixing systems. Size 1000 g, capacity 1260 ml, dimensions 128 x 123 mm (H x outer Ø), jar made in PE, white, with red polypropylene screw lid and anti-tamper ring, bearing aponorm® logo. **Order no. 035263**

**Mixing tool for aponorm® jars for stock preparation
(TOPITEC® jars for stock preparation) SU 1, order no. 025702**

The mixing tool is made in stainless steel. Its shaft is protected by an extra resistant TIN alloy coat. The mixing element (Ø x H = 113.5 x 4.0 mm) is equipped with a wear-proof plastic (PTFE) coat.

**Mixing lid for aponorm® jar for stock preparation
(TOPITEC® jar for stock preparation)**

The mixing lid is made of white POM, with a central tool guide and two recessed O-ring seals. SU = 2, **order no. 025705 (white)**

Replacement seals for mixing lid

SU 6, **order no. 025683**

The silicone seals (Ø x height = 8.1 x 1.6 mm) must always be placed in pairs in the central opening of the mixing lid.

Tool shaft

with titanium-nitride (TiN) alloy coating, in storage box

order no. 025680

TOPITEC® jar holder (size 1)

for TOPITEC® jar sizes 20 g and 30 g

order no. 026631

TOPITEC® jar holder (size 2)

for TOPITEC® jar sizes 50 g to 200 g

order no. 026632

Single-patient jar (complete)

consisting of white production jar and white filling plunger made of POM, with a threaded M9 adapter made of stainless steel screwed in for tube threads, can be sterilised. With storage box (order no. 085500) and user instructions (order no. 080897)

Size 50 g: **order no. 025707**

Size 100 g: **order no. 025696**

Size 200 g: **order no. 025669**

Replacement lift bases

made in white PE, SU=10

order no. 025668

9. MAINTENANCE / CLEANING / INSPECTION INTERVALS & FUNCTION TESTING

Warning

Chapter "**Maintenance / cleaning / inspection intervals & function testing**" contains important information for

- maintenance and repair work, and tests to be performed by electricians at the prescribed inspection intervals.
- Cleaning work might also be carried out by instructed personnel.

Electrician

A person who has the necessary professional training and experience to identify and prevent risks.

The TOPITEC® AUTOMATIC has been designed specifically for use in pharmacies and is manufactured at our own production plants. Our technical expertise and experience enable us to offer you reliable high-tech equipment.

To ensure trouble-free operation over its service life, strictly adhere to our instructions for cleaning, maintenance and operation.

Warning

Operating and maintenance personnel are trained at the site where the TOPITEC® AUTOMATIC is installed. If you have any queries, please contact your dealer.

It is essential that the system is cleaned and maintained at regular intervals in order to enable fault free operation of the TOPITEC® AUTOMATIC mixing system.

During operation, the TOPITEC® AUTOMATIC is exposed to vibration so that screws and clamps might become loose over time. To prevent damage, regularly check the TOPITEC® AUTOMATIC for loose connections.

Protect the device against mechanical impact!

Do not drop the device or any of its accessories. When moving the device on the worktop, ensure that no electronic or mechanical components are damaged.

Protect the device against dust!

Install and store the system and its accessories in a dust-free environment, as fine particles could damage moving parts!

Protect the device against high temperature!

Do not install or store the system or its accessories in a room where they are exposed to high temperatures. High temperature could damage electronic components, or lead to the deformation or melting of certain plastic parts.

Protect the device against extreme temperature changes!

Do not install or store the system or its accessories in a room where they are exposed to very low temperatures. As, during operation, the device is heated to operating temperature, there is a risk of condensation inside the housing, which could damage electronic components.

Note

The average service life of the TOPITEC® AUTOMATIC is 10 years. After this time has elapsed, the device must undergo a complete service by a specialist workshop.

Proper shut-down procedure:

Danger

Before carrying out any cleaning, maintenance or repair task, shut down the device, following the **proper shut-down procedure** (see chapter 4.5)!

Cleaning of housing and base

Do not use pointed implements or aggressive detergents. Clean the housing and the device base with a moist, soft cloth and a mild soapy solution. Alternatively, use an alcohol/water mixture.

To prevent residue, we recommend wiping the device dry with a soft cloth.

Before cleaning the device, disconnect the power plug from the socket!

Cleaning of TOPITEC® AUTOMATIC tool shaft

Step 1: After use, wipe the shaft with a dry cellulose cloth/tissue paper (dry cleaning).

Step 2: Use a cellulose cloth which has already been soaked with a surface disinfectant (e.g. WEPA surface disinfectant 1,000 ml with spray head, order no. 040141) or 2 propanol 70% for thorough cleaning. You can of course also use disinfecting wet wipes such as Meliseptol® HBV. Dispose of the used wipes/tissue paper.

Step 3: Store the tool shaft in the storage box supplied.

The tool shaft can also be cleaned in a rinsing bath, provided that only its bottom part (mixing disc holder) is immersed in water! After cleaning, dry the shaft.

You can find further information at:

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topitec.de/Support/Support-Video

Statutory requirements re. cleaning

Please draw up internal cleaning and hygiene regulations that take into account the relevant statutory requirements for pharmacies, e.g. ApBetrO, BAK Guidelines, GD Hygiene Guidelines, PSA, trade regulations, etc.

9.1 Maintenance

9.1.1 General maintenance instructions

- Correct maintenance at the prescribed intervals is a prerequisite for trouble-free operation of the mixing system. The maintenance intervals in this manual are based on information provided by the manufacturers of third-party components and the experience of WEPA Apothekenbedarf GmbH & Co KG.

9.1.2 Inspection

- **Before starting the device, check the power cable and the main switch for proper connection/functioning** (insulation free of defects, no visible wires).

9.2 Cleaning

Danger

Before carrying out any cleaning, maintenance or repair task, **shut down the device**, following the proper shut-down procedure (see chapter 4.5). Clean the device and the accessories only with implements and tools that have been specifically devised for this task.

9.3 Inspection intervals-function testing

In order to guarantee seamless and safe operation of your TOPITEC® AUTOMATIC mixing system, it is essential that you check your mixing system for external damage at regular intervals and also check switches and buttons for proper functioning. With one-shift operation, perform a function test of all switches and buttons every three months. Make sure that the warning signs on the device are visible at all times. Before starting the device, check the power cable and the main switch for proper connection/functioning. This check must be repeated at least once a year. All rotating and moving device components must be inspected at least once a month. If you encounter any problems or have any queries, please contact your dealer.

9.4 Warning signs attached to TOPITEC® AUTOMATIC mixing system

The following warning signs are attached to the TOPITEC® AUTOMATIC when delivered; Check these signs regularly to ensure that they are properly visible:

Warning	Position on device
	<ul style="list-style-type: none">• Power connection
	<ul style="list-style-type: none">• Base plate
	<ul style="list-style-type: none">• Base plate

Table 3 Warning signs attached to TOPITEC® AUTOMATIC

9.5 Warranty and liability

WEPA Apothekenbedarf GmbH & Co KG (hereafter referred to as "the Manufacturer") grants the end customer (hereafter referred to as "the Customer") based in ... a 3-year warranty from the date of delivery to the customer for the TOPITEC® AUTOMATIC mixing device (subsequently referred to as "Product") stating that the product will be free of material or processing faults for this period (warranty period). The Manufacturer reserves the right to choose between a repair and the supply of new or fully reconditioned parts in relation to faults alleged in connection with this warranty. The warranty terms do not in any way affect the Customer's contractual and statutory rights.

Not covered by warranty are:

- Damage or wear to the Product caused by improper or incorrect operation that is in contravention of best practice and the instructions of the Manufacturer (as laid down in this operating manual)
- Products that show signs of interference or (attempted) repair by a workshop that is not authorised by the Manufacturer to carry out such repairs
- Products that have been modified or operated with unapproved accessories
- Products whose serial number has been removed or made illegible on the type plate

To make a claim under warranty, the Customer must notify the Manufacturer prior to returning the Product by calling **+49 (0)2624 107-361** or writing to **info@wepa-apothekenbedarf.de**. The Customer must allow the Manufacturer a period of 5 working days for the performance of a preliminary fault analysis (by telephone). Claims under this warranty shall only be considered after the Product has been returned to the Manufacturer to the following address:

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Customers based outside Germany may return the Product to their local dealer. To make a claim under warranty, please also furnish us with the original invoice bearing the date of purchase. The costs for the return of the Product shall be covered by the Manufacturer.

If a claim is made under warranty, and the inspection of the Product by the Manufacturer reveals that there is no fault, or that the damage is not covered by warranty for one of the reasons listed above, the Manufacturer shall be entitled to charge a reasonable service fee. This does not apply in cases where the Customer can prove that he could not be aware of the fact that the warranty had been voided.

The Manufacturer warranty as outlined above, including all terms and conditions, shall be transferrable to any new owner based in ..., provided that the original purchase invoice is submitted.

This warranty is governed by German law.

10 PRACTICAL TIPS – INSTANT REMEDY!

When the STOP button is pressed,

message "Stop: program abort" is displayed on the display. Press the **(Stop)** button and switch off the TOPITEC® AUTOMATIC at the power switch. Eliminate the cause of the problem.

Switch on the system.

The display shows two alternating messages: "Release tool shaft" and "Move slide". After having released the shaft locked and moved the slide by hand, the TOPITEC® AUTOMATIC is again in its initial position and ready for operation. The display shows the default settings (M: to 200 g; 0:00 Min; 000 rpm).

Select the desired mode (mixing, grinding, powder-mixing)
Or select the required "Prep. No." using the **(+)** or **(-)** button.
You can now continue working with the TOPITEC® AUTOMATIC system.

Locking device is released during production

Repeat opening and closing of the two locking handles of the clamping device might lead to a reduction of the clamping force. In this case, the locking device might begin to move up during production.

The display texts guide you step by step through the process.

DISPLAY TEXTS:

"Safety bracket tripped!"

REMEDY:

Reposition the safety bracket and follow the instructions on the display.

"Lifting motor monitoring"

Switch off and restart. Please contact your sales partner if the fault continues to occur.

"Torque motor monitoring!"

Switch off and restart. Please contact your sales partner if the fault continues to occur.

"Limit switch error. program abort!"

Switch off and restart. Please contact your sales partner if the fault continues to occur.

11. EMERGENCY-STOP

Operate the power switch and disconnect the power cable in an emergency.

12. DISPOSAL

Disposal

The system, except for the electronic equipment, is mainly made from steel, with some parts in aluminium. Dispose of it through recycling **according to the** applicable statutory waste disposal regulations.

These regulations are likely to require the separate handling and disposal of

- electronic and electrical waste (printed circuit boards);
- batteries, fluorescent lamps/energy-saving bulbs, hazardous substances;
- plastics (housing);
- sheet metal, steel, copper, aluminium (separated by metal).

Before disposing of the device, decontaminate all parts that have been in contact with process substances. Remove any hazardous material from the mixing system.

All oils, solutions, cleaning agents and contaminated cleaning tools (e.g. brushes, cloths) must be disposed of in keeping with the local requirements in accordance with the applicable waste disposal regulations, and in compliance with the information in the material safety data sheets of the manufacturer.

Batteries:

Do not dispose of the batteries through normal household waste. You are legally obliged to ensure that batteries are processed by a specialist disposal company. For this purpose, you can take your old batteries to the public collection points in your district free of charge or anywhere where batteries are sold. Alternatively, you can send the batteries to us (for address, see page 3). We will then dispose of them properly and free of charge for you. Batteries containing toxic substances are marked with a symbol, consisting of a crossed out dustbin and the chemical symbol (Cd, Hg or Pb) of the heavy metal crucial for the toxic classification. The crossed-out waste bin symbol indicates the following: Batteries must not be discarded in household waste. The chemical symbols indicate the following: Pb: battery contains lead; Cd: battery contains cadmium; , Hg: Battery contains mercury.

14. INDEX

Connection of peripherals	31
aponorm® jars / TOPITEC® jars	48-50
Switching off	34
Control elements	33
Operation	30-47
Operating personnel	10
Printer connection port	30-31
Switching on	34
Electrician	18
Disposal	62
Commissioning	26-29
Warranty	59
Safety warnings	15-16
Noise emission	13
Basic equipment	26
Troubleshooting	60-63
Content	4
Emergency-stop	62
Duties of device owner	19-20
Practical tips	60-61
Cleaning	19, 22, 25, 30, 34, 54, 56
Directives and guidelines	2, 19
Interfaces	30
Service	8
Safety bracket	32
Safety precautions	18-20

Power supply	13
Pricing	48-49
Technical data	12
TOPITEC® jars / aponorm® jars	48-50
Type plate	16
Diagram of TOPITEC® AUTOMATIC	9
Maintenance	54, 55, 57, 58
Accessories	50-52

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4. Auflage: Dezember 2020 © Copyright 2013
4th edition: December 2020 © Copyright 2013
4ème édition : décembre 2020 © Copyright 2013
4a edizione: dicembre 2020 © Copyright 2013
4e editie: december 2020 © Copyright 2013
4 naklada: December 2020 © Copyright 2013